DNBSEQ[™] SERVICE OVERVIEW ChIP Sequencing

BG

We care for your samples from the start through

to the result reporting. Highly experienced

laboratory professionals follow strict quality

procedures to ensure high quality results.

Project Workflow

Service Description

ChIP Sequencing is widely used to analyze protein interactions with DNA. It combines chromatin immunoprecipitation (ChIP) with massively parallel DNA sequencing to identify binding sites of DNA-associated proteins, and can be used to precisely map global binding sites for any protein of interest. ChIP sequencing offers higher resolution and more precise and abundant information in comparison with array-based ChIP-chip. Besides clean sequencing data output, BGI offers standard, advanced and customized bioinformatics services to suit your specific research needs.

Confidence:Great correlation with HiSeq data.Low input:As low as 5ng ChIP-ed DNA/sample for human sample.Comprehensive analysis:Correlation analysis between ChIP-Seq and RNA-Seq.

Sequencing Service Specification

The BGI ChIP-Seq Service will be performed with BGI's DNBSEQ sequencing technology, featuring combinatorial probe-anchor synthesis (cPAS) and DNA Nanoballs (DNB) technolog^(j) for superior data quality.



Services

- 50bp Single-end sequencing reads
- Standard output 20 Million reads per sample
- Clean data and bioinformatics analysis are available in standard file formats
- · Available data storage and bioinformatics applications
- · Cloud-based data storage and delivery system



Quality Data

Turn Around Time

- Typical 25 working days from sample QC acceptance to final data delivery
- Expedited services are available, contact your local BGI specialist for details

€

Fast TAT

Cost Effective



Data Analysis

In addition to clean data output, BGI offers a range of standard, advanced and customized bioinformatics pipelines for your ChIP-Seq analysis project, including the correlation analysis of differential expression genes and peak-related genes. Reports and output data files are delivered in industry standard FASTQ, and Excel file formats with publication-ready tables and figures.

STANDARD BIOINFORMATICS ANALYSIS

- Data filtering
- Alignment to the reference genome
- Peak scanning and annotation
- · Identification of differential peaks between samples
- Identification of differential peaks between groups
- · Differential peaks annotation

ADVANCED BIOINFORMATICS ANALYSIS

Motif Analysis

CUSTOMIZED BIOINFORMATICS ANALYSIS

- · Statistics of the epigenetic modification level and mRNA expression level of related genes
- The distribution of epigenetic modification level in different gene categories
- Overall connection of epigenetic modification level and the mRNA expression level in different gene categories
- The relationship of the ratio of epigenetic modification and the ratio of mRNA expression in a pair of samples
- · Clustering analysis based on epigenetic modification level and mRNA expression level
- · Calculate different level of epigenetic modification when mRNA expression level is different
- GO, Pathway analysis, related functional excavation and verification of the genes that differences exist in both epigenetic modification and mRNA expression
- Further customization of Bioinformatics analysis to suit your unique project is available:
- · Please contact your BGI technical representative.

DNBSEQ sequencing technology

DNBSEQ is an innovative high-throughput sequencing solution, developed by BGI's Complete Genomics subsidiary in Silicon Valley. The system is powered by combinatorial Probe-Anchor Synthesis (cPAS), linear isothermal Rolling-Circle Replication and DNA Nanoballs (DNB[™]) technology, followed by high-resolution digital imaging.

The combination of linear amplification and DNB technology reduces the error rate while enhancing the signal. The size of the DNB is controlled in such a way that only one DNB is bound per active site on the DNBSEQ flow cell. This densely patterned array technology provides optimal sequencing accuracy and increases flow cell utilization.



Sequencing data performance

Excellent agreement of data quality between HiSeq and DNBSEQ

The plot below shows the high sequencing data quality of human HeLa cell libraries and mouse libraries on DNBSEQ as compared to the HiSeq platform. By applying the same filtering criteria, the clean rate and unique mapping rate from the same samples are consistent on both platforms.



Excellent agreement of reads per million (RPM) between HiSeq and DNBSEQ

The scatterplots below compare the reads per million (RPM) values of human HeLa cell sample sequenced on the DNBSEQ and HiSeq platforms. The coefficients of sequencing repeatability on both platforms are high (Figures a, b). There is a high correlation between DNBseq and HiSeq platforms as well (Figure c).



Excellent agreement of peak number detection between HiSeq and DNBSEQ

The figures below show the comparison of detected peak numbers on both HiSeq and DNBSEQ platforms from the same samples^[2]. The common peak detection rate of both drosophila and human HeLa cell sample are higher than 99%, while both platforms have a small amount of their own uniquely detected peaks.



Sample Requirements

We can process DNA samples of human, plant, animal and microbial samples, with the following general requirements:

	DNA Amount and Concentration	Minimum Volume
ChIP-ed DNA	Amount \ge 10ng, Concentration \ge 1 ng/µl	15 µl
ChIP-ed DNA (human sample only)	Amount \ge 5ng, Concentration \ge 1 ng/µl	15 µl

References

[1] Human genome sequencing using unchained base reads on self-assembling DNA nanoarrays
Drmanac R, Sparks AB, Callow MJ, Halpern AL, Burns NL, Kermani BG, Carnevali P, Nazarenko I, Nilsen GB, Yeung G, et al.
Science (2010) 327(5961):78–81 doi:10.1126/science.1181498
http://science.sciencemag.org/content/327/5961/78.full
[2] ChIP-seq guidelines and practices of the ENCODE and modENCODE consortia
Stephen G. Landt,Georgi K. Marinov,Anshul Kundaje, et al.
Genome Res. 2012 Sep;22(9):1813-31. doi: 10.1101/gr.136184.111
https://www.ncbi.nlm.nih.gov/pubmed/22955991



Want to learn more?

Contact your local BGI account representative for the most affordable ChIP-sequencing rates in the industry, to discuss how we can meet your specific project requirements or for expert advice on experiment design, from sample to bioinformatics.

info@bgi.com www.bgi.com

BGI Offices

BGI Americas

One Broadway, 14th Floor Cambridge, MA 02142, USA

BGI Europe Ole Maaløes Vej 3, DK-2200 Copenhagen N, Denmark

BGI Asia

Building NO.7, BGI Park, No.21 Hongan 3rd Street, Yantian District, Shenzhen

For Research Use Only. Not for use in diagnostic procedures (except as specifically noted). Copyright© BGI 2022. All trademarks are the property of BGI, or their respective owners. This material contains information on products which is targeted to a wide range of audiences and could contain product details or information otherwise not accessible or valid in your country. Please be aware that we do not take any responsibility for accessing such information which may not comply with any legal process, regulation, registration or usage in the country of your origin. Note, BGI's genetic testing products have not been cleared or approved by the US FDA and are not available in the USA. For Research Use Only. Unless otherwise informed, all sequencers and sequencing reagents are not available in Germany, USA, Spain, UK, Hong Kong, Sweden, Belgium or Italy. Certain sequencing services are not available in USA and Hong Kong. Please contact a representative for regional availability. The company reserves the right of final interpretation.

BGI Australia

L6, CBCRC, 300 Herston Road, Herston, Brisbane, Queensland 4006, Australia



